

Rotary Mill

for grinding a variety of materials for subsequent analyses



Brabender®

... where quality is measured.

Brabender[®] Rotary Mill

for grinding a wide variety of materials for subsequent laboratory analyses


The material is fed into the rotary mill's milling chamber via a feed funnel. It is then dosed via a slider on the funnel's outlet.

The mill's interior features four fixed blades made from a special steel. The rotor fixed to the motor shaft is fitted with six interchangeable, adjustable blades, whose cutting edges work against those of the fixed blades, shredding the material. The shredded material then falls through the sieve into the collection container.

All residues are milled further, until the desired degree of fineness has been achieved.

Technical Data

- Dimensions (BxHxT): 320 x 600 x 610 mm
- Weight: ca. 55 kg
- Mains connection: 3x 400 V; 50/60 Hz + N +PE;
2,6 A 3x 230 V; 50/60 Hz + PE; 4,7 A

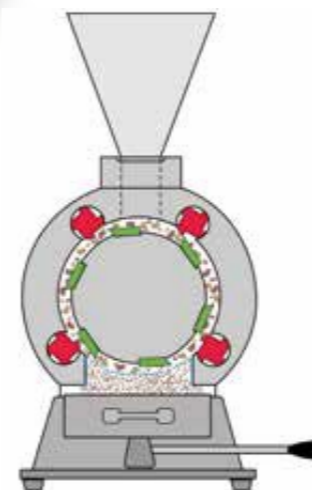

Rotary Mill (schematic diagram)

Highlights

Knives made of high-quality special steel

- 4 fixed knives in the grinding chamber, 6 exchangeable and adjustable knives on the rotor (see schematic diagram)

Interchangeable sieves

- with different finenesses (perforations from 0.5-5.0 mm available)

Low cleaning effort

- Dust leakage is significantly reduced thanks to an effective sealing system

Safety provisions

- Microswitch on the door, safety switch on the collecting hopper and in the grinding housing

Application examples

- Grains (analyses according to Farinograph whole grain method (ICC standard no. 15/1), Falling Number, Glutomatic)
- Plastics
- Tobacco
- Fibrous materials (hay, grass, straw, leaves etc.)
- Tough materials (leather, animal skins, linoleum, pulp etc.)
- Solid materials (pasta, roots, coconut shells etc.)

Brabender® GmbH & Co. KG

Headquarters Germany
Kulturstrasse 49 - 51
D-47055 Duisburg
Phone: +49 203 7788 0
sales@brabender.com
www.brabender.com

C.W. Brabender® Instruments, Inc.

USA
50 East Wesley Street
South Hackensack, NJ 07606
Phone: +1 201 343 8425
sales@cwbrabender.com
www.cwbrabender.com

000 Brabender®

Russia
Yagodinskaya St. 25
420032 Kazan
Phone: +7 843 233 46 66
000.brabender@brabender.ru
www.brabender.ru



Headquarter Germany



Contact us 24/7: +49 203 7788-131

Brabender agencies worldwide.
© 2021 Brabender GmbH & Co. KG

All trademarks are registered. Subject to design
and technical modifications without notice.

Brabender®
... where quality is measured.